

## I200 Two-channel Digital Electrometer with servo control option

### Features

- Two independent gated integrator channels
- Dynamic range 0.1pA to 100uA
- Integrated digitization and communications
- Integrated calibration test source
- Selection of current and charge integration modes
- External trigger capability
- Optional high-precision analog output for servo control applications
- Auxiliary I/O port

### Applications

- Ionization chambers
- Two electrode beam position monitors
- Photodiodes
- Ultra-low current and charge measurement
- With servo control option:
  - Stabilisation of double crystal monochromators for synchrotron beamline experiments by control of the crystal cage piezo motor amplifier
  - Stabilisation of beam position in charged particle or photon beamlines



### Options

- Auxiliary HV output 0 to 1000V
- Alternative feedback capacitor options
- Servo control options

### Specifications


Operating principle	Gated integrator (charge integrating amplifier)
Integration capacitor	Dual, selectable
Input noise	< 10 fA rms + 1 fA rms per pF input load up to 100 pF (1 second integration, 10 pF capacitor, 0 V bias)
Input offset current	< 100 fA at 0V bias, 25 C < 600 fA at 0V bias, 50 C < 200 fA at 400V bias, 25 C
Stability Output drift	< 5 ppm / C / hour
Integration time	User selectable, 100 µsec minimum
Integration modes	0: continuous current; 1: continuous charge, 2: lossless charge; 3: external gate integration; 4: externally-triggered burst mode
External gate	0/+5 V, 50 ohm or fiber-optic receiver, software-selectable
Digitization	16 bit bipolar, 250kHz
Auxiliary HV PSU	(Factory option) 0 to 1000V programmable (polarity and maximum voltage factory selectable), 1mA max. Noise and ripple < 0.1%
Power input	+24V (+/- 2V) DC, 350mA typ, 500mA max
Controls	Two rotary switches for loop address and comms mode/baud rate
Displays	Status LEDs (power, device status, comms mode, data transmission rcv/xmit). "HV on" LED
Case material	Stainless steel sheet
Weight	1.64kg (3.6 lb)
Operating environment	10 to 35C (15 to 25 C recommended to reduce drift and offset) , < 70% humidity, non-condensing, vibration < 0.1g all axes (1 to 1000Hz)
Shipping and storage environment	-10 to 50C, < 80% humidity, non-condensing, vibration < 2g all axes, 1 to 1000Hz Hz

## Interfacing

Interfaces	RS-232, 8-bit ASCII. Selectable baud rate USB, 8-bit ASCII 3 Mbit/sec Fiber-optic loop, 10 Mbit/sec serial, 9-bit asynchronous binary. Ethernet connection to host through A300 or A500 loop controllers
Host computer	ASCII communications based on SCPI. Diagnostic host program supplied for Microsoft® .net framework. DLLs available for Microsoft® .net, National Instruments™ Labview™ and Microsoft® C++

## Servo control option

Principle	The I200 measures one or two currents, from devices such as ionisation chambers, photodiodes, or secondary electron emission electrodes. The value of an arithmetic function of the measured currents is maintained by adjusting a 0 to 10 V control output using a PID control algorithm
Typical applications	Beam intensity stabilization for Double Crystal Monochromators Beam position control in charged particle or photon beamlines


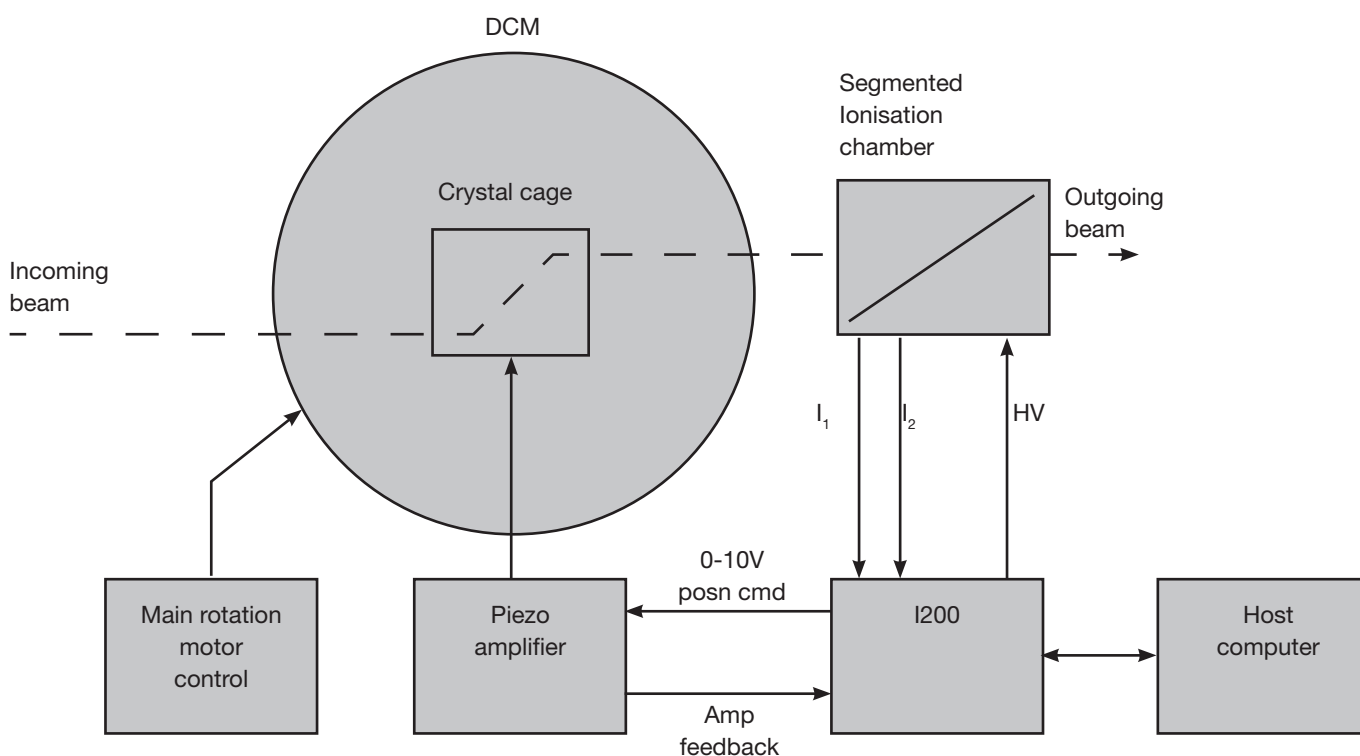


## Sample application - double crystal monochromator stabilisation


Servo type	PID controller implemented in firmware; servo parameters software configurable
Process variable choices	$I_1$ , $(I_1 / I_2)$ , $(I_1 + I_2)$ , $(I_1 - I_2)$ , $((I_1 - I_2) / (I_1 + I_2))$
Analog output	16 bit DAC, 0 to +10 VDC
Analog input	10 bit ADC, 0 to +10 VDC
Servo frequency	200 Hz maximum
Automated functions	Open loop direct control of the DAC output Peak scan Peak spectrum upload Servo suspension if process variable bad

# Connectors

Signal inputs	Two triaxial three lug bayonet. Mates with Trompeter PL74-7 or equivalent
Auxiliary HV out	SHV
External gate in	a) BNC (isolated from case) b) Avago HFBR ST bayonet
Auxiliary I/O and servo control	9 pin D female 1 Analog Gnd      6 Analog output 2 Analog Gnd      7 Analog input 3 +24 VDC out    8 PSU Gnd 4 Opto common    9 Opto in B 5 Opto in A
USB	USB B type female
RS-232	Six pin mini-DIN ("PS/2") 1 Tx                    4 n/c 2 Rx                    5 n/c 3 Gnd                  6 n/c
Fibre optics	Two 1mm Avago HFBR ST bayonet
Power in	2.1mm threaded jack. Mates with Switchcraft S761K or equivalent
Ground	M4 threaded stud

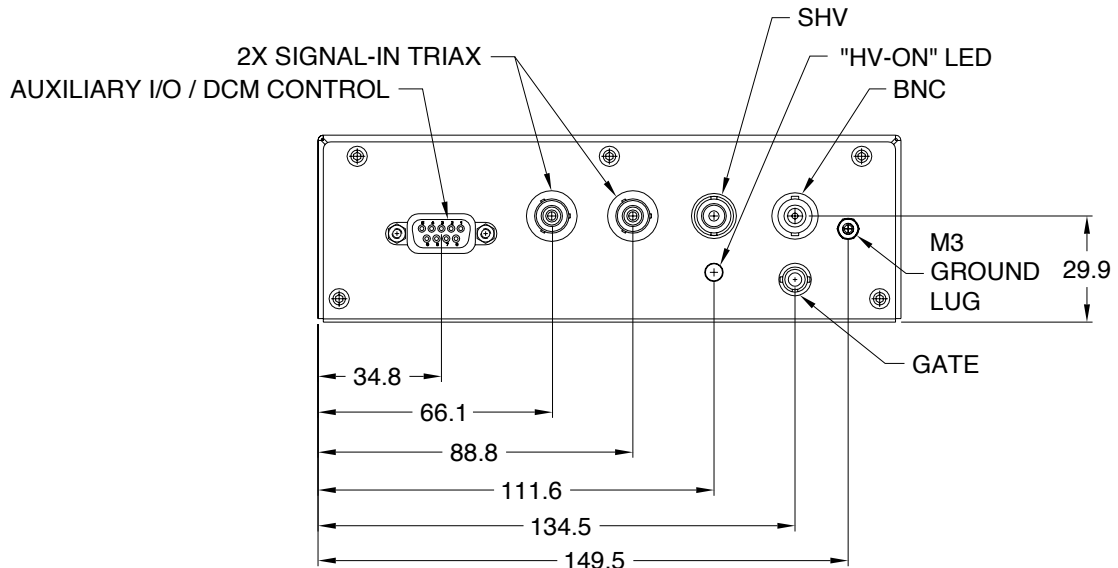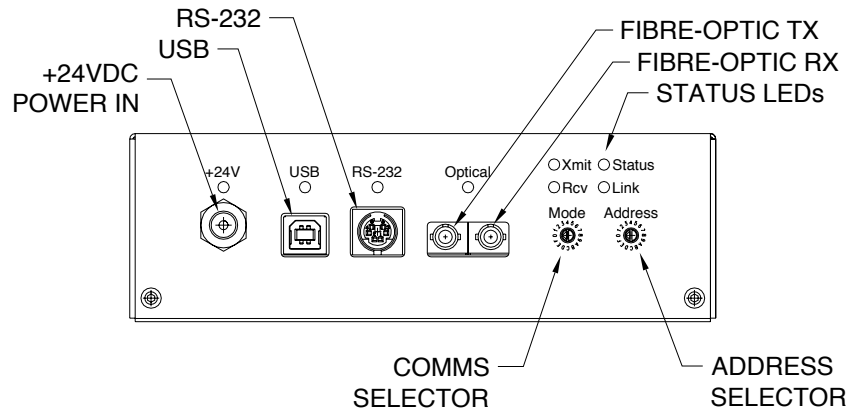

Dims mm

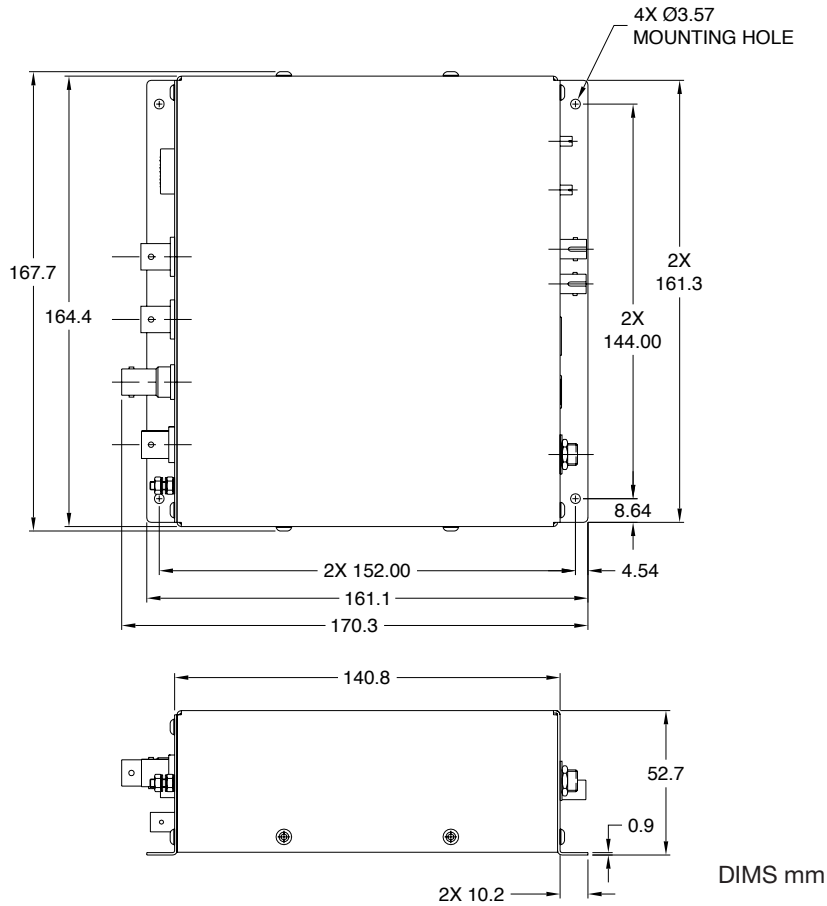

## Ordering information

I200 two channel electrometer with user manuals, software drivers, calibration data

Available on request

Add auxiliary HV supply positive 1000 V / 500 V / 200V

Specify channel 1 feedback capacitors x pF, y pF (default is 10pF, 1000pF)

Specify channel 2 feedback capacitors x pF, y pF (default is 10pF, 1000pF)

Add multimode servo control option

Add modulation mode servo option

Your local agent is:

**Head Office**  
 FMB Oxford  
 Unit 1 Ferry Mills  
 Osney Mead  
 Oxford OX2 0ES  
 United Kingdom

Tel +44 (0)1865 320300  
 Fax +44 (0)1865 320301

[enquiries@fmb-oxford.com](mailto:enquiries@fmb-oxford.com)



[www.fmb-oxford.com](http://www.fmb-oxford.com)